

FFPE Tissue Genomic DNA Extraction Kit

Instruction for Use (V1.1)

[REF] UBE-S00450Q

[Specification] 50 preps/kit

[Research Use]

Ultrassay FFPE Tissue Genomic DNA Kit is optimized to extract genomic DNA from formalin-fixed and paraffin-embedded tissues. The kit uses specific lysis buffer to release DNA from FFPE tissues. DNA binds to silica-based column and contaminants flow through. DNA is eluted from the membrane. The isolated DNA is suitable for PCR, qPCR and other molecular biology applications.

- Simple and fast
- Complete remove of contaminants and inhibitors

[Kit Contents]

No.	Content	50 preps
1	Lysis Buffer 15 (LB15)	11ml
2	Binding Buffer 15 (BB15)	11ml
3	Clean Buffer 15 (CB15)	6 ml
4	Wash Buffer 15 (WB15)	12 ml
5	Elution Buffer (EB)	25 ml
6	Proteinase K (20 mg/ml)	1 ml
7	Genomic Spin Column with Collection Tubes	50 pcs

[Storage Conditions and Shelf Life]

At room temperature (15-25°C) for up to 12 months.

[Application]

PCR, restriction enzyme digestion, and other downstream applications.

[Test Procedures]

Before use, add 100% ethanol to CB15 (24ml) and WB15 (48ml). Prepare 56°C water bath or heating block.

(All centrifugation steps are performed at room temperature)

1. Preparing Materials

A. Formalin-fixed tissue: Place 10-30 mg chopped tissue into a sterile 1.5 ml microcentrifuge tube.

- Add 500 µl 1×PBS and mix thoroughly by vortex. Incubate at room temperature for 1 minute. Centrifuge at 12,000×g for 30 seconds and remove the supernatant.
- Repeat the above step twice.

B. FFPE sample block: Trim excess paraffin off the tissue and scrap 10-30 mg tissue by a scalpel.

FFPE section: Scrap tissue form 3-10 sections (5-10 µm thick) and place into a sterile 1.5 ml microcentrifuge tube.

- Add 1 ml xylene to the sample, close the lid and vortex vigorously for 10 seconds. Centrifuge at 12,000×g for 2 min, then remove the supernatant by pipetting.

(It is recommended to carefully perform this step in a fume hood because xylene is high toxic chemical).

- Add 1 ml ethanol (96–100%) to the pellet, and mix by vortex. Centrifuge at 12,000×g for 2 minutes, then remove the supernatant by pipetting.
- Open the tube and incubate at room temperature or up to 37°C until all residual ethanol has evaporated.

- Add 200 µl of LB15 and 20 µl of Proteinase K to resuspend the pellet, mix thoroughly by vortex. Incubate at 56°C for 1 hour (or until the sample has been completely lysed).

In the meantime, prepare 90°C water bath or heating block for the next step. If the water bath or heating block from this step will be used for the next step, we suggest to leave the sample at room temperature after this incubation and wait until the water bath or heat block has reached 90°C.

- Incubate at 90°C for 1 hour. Briefly centrifuge the tube to remove drops from the lid.

The incubation at 90°C in LB15 Buffer can partially reverse formalin-induced crosslinking of nucleic acids. Longer incubation time and higher temperature may result in more fragmented DNA.

Optional: If RNA-free genomic DNA is needed, add 10 µl of RNase A (20 mg/ml) to the lysate, incubate at room temperature for 2 minutes.

- Add 200 µl of BB15 to the sample, and mix thoroughly by vortex (white precipitates may form).

- Add 250 µl of 100% ethanol, and mix thoroughly by vortex (if there are white precipitates, vortex until they disappear). Briefly centrifuge the 1.5 ml tube and transfer all the lysate to a Genomic Spin Column, then centrifuge the tube at 12,000×g for 1 minute, discard the flow-through.

- Add 500 µl of CB15 (make sure that ethanol has been added), and centrifuge at 12,000×g for 30 seconds, discard the flow-through.
- Add 500 µl of WB15 (make sure that ethanol has been added), and centrifuge at 12,000×g for 30 seconds, discard the flow-through.

- Repeat step 7 once.










- Centrifuge the Spin Column at 15,000 ×g for 2 minutes. Open the tube and air-dry the Spin Column at room temperature to volatilize the residual ethanol.

- Place the Spin Column in a sterile 1.5 ml microcentrifuge tube. Add 30-100 µl of Elution Buffer (for higher yield, prewarm the buffer to 65°C) or distilled water (pH >7.0) to the center of the membrane. Incubate at room temperature for 1 minute. Centrifuge at 12,000×g for 1 minute to elute the DNA.

Notes

- Too many starting materials will affect DNA yield.
- Fix tissue samples as quickly as possible after surgical removal. Use a fixation time of 14-24 hours (longer fixation times lead to more fragmented DNA). Thoroughly dehydrate samples prior to embedding because residual formalin can inhibit the Proteinase K digestion.
- Use sterile tubes and pipette tips to avoid DNase contamination.

[Index of Symbols]

Symbols	Meanings	Symbols	Meanings
	Use By		Date of Manufacture
	Temperature Limitation		Consult Instructions for Use
	Lot Number		Manufacturer
	Number of Tests		Reference Number
			Any warnings and/or precaution to take



Ultrassay Biotech Co., Ltd.

Add: 2906, Building B, Dongyi Square, No. 169 Funan Road, Hefei, Anhui, China

Tel: +86-551-6288 1663

Website: [Http://ultrassay.com](http://ultrassay.com)

Version: A/0 Issue date: 2021.12.21